







DEOXYRIBONUCLEASE I

Molecular Biology/Cell Biology

Worthington prepares Deoxyribonuclease I from bovine pancreas at different levels of purity to suit many different applications. Product Codes: DPRF and DPRFS are especially designed for Molecular Biology applications and contain the lowest levels of ribonuclease and protease activity. They are suitable for use in techniques requiring digestion of DNA in the recovery of intact RNA or where the integrity of structural proteins or enzymes must be maintained. Applications have included nick translation, DNA mapping, isolation of nuclear RNA and protein, plasmid construction, and RNA polymerase synthesis of RNA. Additional grades for other applications are listed below.

Recombinant DNase I, Animal Free (AF)

Bovine pancreas is a rich source of RNase which is often found in many commercial DNase preparations. Producing DNase I by recombinant means in an organism with much lower levels of endogenous RNase greatly facilitates purification of an enzyme with undetectable levels of RNase (DR1/DR1S). The processes involved in the production and isolation of recombinant DNase I are completely devoid of animal based components which eliminates the possibility of introducing animal derived pathogens into bioprocessing procedures. Recombinant DNase I is suitable for such applications as:

- Removing genomic DNA from RNA preparations prior to RT-PCR
- Degradation of DNA templates after transcription reactions
- Removing unwanted DNA from samples prior to Northern blotting
- Removing DNA during biopharma and bioprocessing procedures

Produced in Pichia pastoris, Ribonuclease & Protease Free, Lyophilized Powder.Units per mg proteinLS006362 LS00636050 ku735.Molecular Biology Grade. Free of RNase and protease. Chromatographically purified and lyophilized powder containing glycine as a stabilizer. Store at 2 - 8°C.>2 Kunitz units per microliterDR1SLS006353 LS0063552 ku\$ 47.DNase I, Recombinant (AF), Produced in Pichia pastoris, Ribonuclease & Protease Free, Solution.>2 Kunitz units per microliterDR1SLS006353 LS0063572 ku\$ 47. 5x2 kuNo I e cularBiology Grade. Chromatographically purified as a ready-to-use solution at ≥ 2 Kunitz U/µl in 5mM calcium acetate, 4mg/ml glycine, pH 5.0 and 50% glycerol. Includes IOX reaction buffer. Store at -20°C. REQUIRES ICE PACK.≥ 2,000 Kunitz units per mg dry weightDR2LS006320 LS006323 LS00632325 ku\$ 85. 240. 240.DNase I, Recombinant (AF), Produced in Pichia pastoris, Bioprocess Grade. Chromatographically purified AF bioprocessing grade. Supplied as a lyophilized powder containing glycine as a stabilizer. For the removal of DNA in bioprocessing and primary stem cell isolation applications. May≥ 2,000 Kunitz units per mg dry weightDR2LS006320 LS006323 LS006323 LS00632525 ku S0 ku S00 ku S00 ku\$ 85. 240. S00 ku	Description	Activity	Code	Cat. No.	Size	Price
Drase I, Recombinant (AF), Produced in Pichia pastoris, Ribonuclease & Protease Free, Solution.Link units per microliterLiso063555x2 ku188.Molecular Biology Grade. Chromatographically purified to remove RNase and protease. Supplied as a ready-to-use solution at ≥ 2 Kunitz u/µl in 5mM calcium acetate, 4mg/ml glycine, pH 5.0 and 50% glycerol. Includes 10X reaction buffer. Store at -20°C. REQUIRES ICE PACK. $\geq 2,000$ Kunitz units per mg dry weightDR2LS006320 LS006322 LS006323 LS006323 S00 ku25 ku \$ 85. LS006323 S00 ku Produced in Pichia pastoris, Bioprocess Grade. Chromatographically purified AF bioprocessing grade. Supplied as a lyophilized powder containing glycine as a stabilizer. For the removal of DNA in bioprocessing and primary stem cell isolation applications. May $\geq 2,000$ Kunitz units per mg dry weightDR2LS006320 LS006323 LS00632525 ku Bulk $\geq 8.$	Produced in Pichia pastoris, Ribonuclease & Protease Free, Lyophilized Powder. Molecular Biology Grade. Free of RNase and protease. Chromatographically purified and lyophilized powder containing glycine as a stabilizer. Store at 2 - 8°C.	units per	DR1	LS006362	50 ku	\$ 174.00 735.00 Inquire
Produced in Pichia pastoris, Bioprocess Grade. units per mg dry weight LS006322 100 ku 240. Bioprocess Grade. units per mg dry weight LS006323 500 ku 780. LS006325 Bulk Inqui powder containing glycine as a stabilizer. For the removal of DNA in bioprocessing and primary stem cell isolation applications. May May Image: Containing co	Produced in Pichia pastoris, Ribonuclease & Protease Free, Solution. Molecular Biology Grade. Chromatographically purified to remove RNase and protease. Supplied as a ready-to-use solution at ≥ 2 Kunitz u/µl in 5mM calcium acetate, 4mg/ml glycine, pH 5.0 and 50% glycerol. Includes 10X reaction buffer. Store at -20°C.	units per	DR1S	LS006355	5x2 ku	\$ 47.00 188.00 Inquire
	Produced in Pichia pastoris, Bioprocess Grade. Chromatographically purified AF bioprocessing grade. Supplied as a lyophilized powder containing glycine as a stabilizer. For the removal of DNA in bioprocessing and	units per	DR2	LS006322 LS006323	100 ku 500 ku	\$ 85.00 240.00 780.00 Inquire (0ver) [01.17]

730 Vassar Avenue, Lakewood, New Jersey 08701 800.445.9603 • 732.942.1660 • Fax: 800.368.3108 • 732.942.9270 Worthington-Biochem.com **ISO9001 Certified**



Description	Activity	Code	Cat. No.	Size	Price				
Deoxyribonuclease I, Ribonuclease & Protease Free, Solution. Purified from bovine pancreas. Molecular Biology Grade. Chromatographically purified to remove RNase and protease. Supplied as a solution at approximately 2 Kunitz units per microliter approximately 1 mg/ml containing 50% glycerol and 1 mM calcium chloride. Store at 2-8°C or -20°C.	≥2,000 Kunitz units per ml	DPRFS	LS006342 LS006344 LS006348	100 un 500 un Bulk	\$ 27.00 70.00 Inquire				
Deoxyribonuclease I, Ribonuclease & Protease Free. Purified from bovine pancreas. Molecular Biology Grade. Chromatographically purified to remove RNase and protease. Lyophilized in vials. Each 10,000 unit vial contains 2 mg glycine, 2 µmoles calcium, and \geq 10,000 units of DNase I. Each 2,500 unit vial contains 0.5 mg glycine, 0.5 µmoles calcium, and \geq 2,500 units of DNase I. Dissolving the entire 10,000 unit vial in 5 ml, or the entire 2,500 unit vial in 1.25 ml, provides the equivalent of a 1 mg/ml solution. (ku = 1000 un). Store at 2-8°C. PROTECT FROM MOISTURE.	≥2,000 Kunitz units per mg dry weight	DPRF	LS006331 LS006333 LS006343 LS006334	2500 un 10 ku 50 ku Bulk	\$ 41.00 148.00 540.00 Inquire				
Deoxyribonuclease I. Purified from bovine pancreas. Chromatographically prepared especially for low (≤0.0005%) RNase A activity content. A lyophilized powder containing glycine as a stabilizer. Store at 2 - 8°C. PROTECT FROM MOISTURE.	≥2,000 Kunitz units per mg dry weight	DPFF	LS006330 LS006328 LS006332	25 ku 125 ku Bulk	\$ 86.00 316.00 Inquire				
Deoxyribonuclease I. Purified from bovine pancreas. Chromatographically purified. A lyophilized powder with glycine as a stabilizer. Store at 2 - 8°C. PROTECT FROM MOISTURE.	≥2,000 Kunitz units per mg dry weight	D	LS002004 LS002006 LS002007 LS002009	5 mg 20 mg 100 mg Bulk	\$ 33.00 76.00 270.00 Inquire				
Deoxyribonuclease I, Filtered. Purified from bovine pancreas. Filtered through a 0.22 micron membrane and lyophilized in vials. Material is not tested for pyrogenicity. Store at 2-8°C. PROTECT FROM MOISTURE.	≥2,000 Kunitz units per mg dry weight	DCLS	LS002058 LS002060	11 mg 25 mg	\$ 92.00 158.00				
Deoxyribonuclease I, Standard Vial. Purified from bovine pancreas. Lyophilized in vials for assay standardization. Labeled to show established activity. Not suitable for assays at neutral pH. Store at 2-8°C.	~2,000 Kunitz units per vial	DSV	LS002173 LS002172	2 ku 5x2 ku	\$ 19.00 53.00				
PDS Kit, DNase Vial. Purified from bovine pancreas. A component of the Papain Dissociation System. This material is 0.22 micron membrane filtered and lyophilized in autoclaved vials. A vial reconstituted with 0.5 ml of EBSS or equivalent yields a solution of 2000 units/ml of deoxyribonuclease (1 mg/ml). Store at 2-8°C.	≥1,000 units per vial	D2	LK003170 LK003172	1 vi 5 vi	\$ 23.00 78.00				
Deoxyribonuclease I. Purified from bovine pancreas. Partially purified. A lyophilized powder. Store at 2-8°C. PROTECT FROM MOISTURE.	≥2,000 Kunitz units per mg dry weight	DP	LS002138 LS002139 LS002140 LS002141	25 mg 100 mg 1 gm Bulk	\$ 37.00 90.00 747.00 Inquire				
Deoxyribonuclease I. Purified from bovine pancreas. Partially purified. A lyophilized powder. Store at 2-8°C. PROTECT FROM MOISTURE.	≥1,250 Kunitz units per mg dry weight	DPB	LS002145 LS002147 LS002149	100 mg 1 gm Bulk	\$ 77.00 550.00 Inquire				
Related Products									
Actin • Albumin, Nuclease-Free • Deoxyribonuclease II • Deoxyribonucleic Acid and Related Products • Histones • Lysozyme • Nuclease, Micrococcal Nuclease, S1 • Phosphatase, Alkaline • Phosphodiesterase I • Phosphodiesterase II • Proteinase K									

Nuclease, S1 • Phosphatase, Alkaline • Phosphodiesterase I • Phosphodiesterase II • Proteinase K Reverse Transcriptase, Recombinant HIV • RNase A • RNase T1 • RNase T2 • Ribonucleic Acid

Complete Catalog, Tissue Dissociation Guide and Enzyme Manual available online at: Worthington-Biochem.com TissueDissociation.com