

Bst DNA Polymerase, Exonuclease Minus 50,000 U/ml



Technical Specifications

<p>Catalog No. 30028-1 10,000 Units (200µl @ 50 U/ µl) Includes 10X DNA Polymerase Buffer B (4 X 1.2 ml)</p>	
<p>Store at –20°C. For Research Use Only. Not for use in Diagnostic Procedures.</p>	

Product Description	Bst DNA Polymerase, Exonuclease Minus, 50,000 units/ml.
Storage Buffer	10 mM Tris-HCl, pH 7.5, 50 mM KCl, 1 mM DTT, 0.1 mM EDTA, 0.1% Triton X-100, and 50% Glycerol.
Stability	Bst DNA Polymerase, Exonuclease Minus is stable for one year from the date received if stored at –20°C.
Recommended Reaction Conditions	50 U Bst DNA Polymerase, Exonuclease Minus; 1X DNA Polymerase Buffer B containing 20 mM Tris-HCl pH 8.8, 10 mM (NH ₄) ₂ SO ₄ , 10 mM KCl, 2 mM MgSO ₄ , and 0.1 % Triton X-100.
Activity Determination	One unit catalyzes the incorporation of 10 nmol of dNTP into acid-insoluble material in 30 minutes at 65°C in 20 mM Tris-HCl pH 8.8 , 10 mM (NH ₄) ₂ SO ₄ , 10 mM KCl, 2 mM MgSO ₄ , 0.1 % Triton X-100, 30 nM M13mp18 ssDNA, 70 nM M13 sequencing primer(-47) 24 mer, 200 µM dGTP, dATP, dTTP, dCTP (a mix of unlabeled and [³³ P]dCTP), and 0.1 mg/ml BSA.
Absence of Endonuclease or Nicking Activity	Incubation of 50 U of Bst DNA Polymerase, Exonuclease Minus with 1 µg of supercoiled pBR322 DNA for 16 hours at 37° and 65°C resulted in no detectable conversion to relaxed or linear forms by agarose gel electrophoresis.
Absence of Exonuclease Activity	Incubation of 50 U of Bst DNA Polymerase, Exonuclease Minus with 1 µg of HindIII-cut lambda DNA for 16 hours at 37° and 65°C resulted in no smearing of bands on agarose gels. Single stranded and double stranded exonuclease activities were tested by incubating 10 µl of enzyme at 50 U/ µl with radiolabeled DNA substrate for one hour at 37° and 65°C, resulting in less than 0.1% release of TCA-soluble counts.
Purity	>99% pure by SDS PAGE. No detectable DNA contamination. 10 µl of enzyme at 50 U/ µl of the sample was tested for <i>E. coli</i> genomic DNA contamination by PCR amplifying with the <i>E. coli</i> 16S ribosomal primers.

Applications

1. DNA sequencing through high GC regions (1, 2)
2. Rapid Sequencing from nanogram amounts of DNA template (3)

Heat Inactivation: 80°C for 20 min.

References

- 1) Griffin, H. and Griffin, A. (1994) *PCR Technology*, 228-229.
- 2) McClary, J. et al. (1991) *J. DNA Sequencing and Mapping*, 1, 173-180.
- 3) Mead, D.A. et al. (1991) *Biotechniques*, 11, 76-87.

Lucigen Corporation 2905 Parmenter St, Middleton, WI 53562 USA
Toll Free: (888) 575-9695 | (608) 831-9011 | FAX: (608) 831-9012
lucigen@lucigen.com www.lucigen.com

Legal & Warranty

Lucigen does not encourage or support the unauthorized or unlicensed use of patented nucleic acid amplification processes. It is the sole responsibility of the buyer to ensure that use of the product does not infringe the patent rights of third parties. If the purchaser is not willing to accept these use limitations, Lucigen Corporation is willing to accept return of the product for a full refund.

Lucigen warrants that its products will conform to the standards stated in its product specification sheets in effect at the time of shipment. Lucigen will replace, free of charge, any product that does not conform to the specifications. This warranty limits Lucigen's liability only to the replacement of the product. THIS WARRANTY IS EXCLUSIVE, AND LUCIGEN MAKES NO OTHER WARRANTY, EXPRESS OR IMPLIED, INCLUDING WITHOUT LIMITATION, ANY IMPLIED WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE. The stated express warranties, and the remedy provided for breach thereof, are in lieu of all other liability or obligations of Lucigen for any damages whatsoever arising out of or in connection with the delivery, use, performance, or the inability to use any of its products. IN NO EVENT SHALL LUCIGEN BE LIABLE UNDER ANY LEGAL THEORY (INCLUDING BUT NOT LIMITED TO CONTRACT, NEGLIGENCE, STRICT LIABILITY IN TORT, OR WARRANTY OF ANY KIND) FOR ANY INDIRECT, SPECIAL, INCIDENTAL, CONSEQUENTIAL, OR EXEMPLARY DAMAGES (INCLUDING BUT NOT LIMITED TO LOST PROFITS). Without limiting the effect of the preceding sentence, Lucigen's maximum liability, if any, shall not exceed the purchase price paid by buyer for the product.